

We supply:

- Living materials
- Culture media
- Frozen materials
- Preserved materials
- Skeletons
- Microscopes
- Wall charts
- Stains and Indicators
- Seeds
- Books
- CD-ROMs

*and lots more in our full
colour catalogue*

Hydra Culture Medium Preparation:

20mls of Southern Biological's Hydra medium concentrate will make 1 litre of useable medium. Make up solution as follows:

Instructions:

- Take 20mls of Hydra Culture Medium concentrate (code CM12) and add 950mls of distilled water.
- Check pH and adjust to 7.8-8.0. Use either dilute NaOH or dilute Acetic Acid. (1% solutions added one drop at a time)
- Make up the volume to 1L with distilled water.

Ref: CM12 20/03/09