

Chlorella Culture Medium

Catalogue No. CM 3

We supply:

- Living materials
- Culture media
- Frozen materials
- Preserved materials
- Skeletons
- Microscopes
- Wall charts
- Stains and Indicators
- Seeds
- Books
- CD-ROMs
- Forensic Equipment

*and lots more in our full
colour catalogue or visit our
website catalogue at
www.southernbiological.com*

Chlorella Culture Medium comprises a sterile chemically defined mineral salts solution. It is desirable to keep any unused medium in the refrigerator at 6-8°C for up to 4 weeks after opening in order to minimize bacterial/fungal contamination which may occur during sub-culturing.

CULTURE MEDIUM

Shake the contents of the bottle to disperse any sediment. Pour 20-25ml of the Chlorella Culture Medium into a sterile 30ml bottle. Add 2-3ml of your live Chlorella culture. Replace the lid on the bottle and maintain at 20-25°C, loosely capped, on a bench which receives bright, but not direct, sunlight. Growth will be far stronger and more rapid in spring/summer than in autumn/winter, and stronger lighting is necessary in the latter seasons. Sub-culture every 6-8 weeks.

Ref: CM 3 25.09.09