

Brine Shrimp Eggs

Catalogue Code: L5.00

We supply:

- Living materials
- Culture media
- Frozen materials
- Preserved materials
- Skeletons
- Microscopes
- Wall charts
- Stains and Indicators
- Seeds
- Books
- CD-ROMs
- Forensic Equipment

*and lots more in our full
colour catalogue*

DIRECTIONS FOR HATCHING:

1. Prepare a salt nutrient solution by adding coarse salt to tap water in a suitable sized container (use 1 teaspoon of salt to 1 teaspoon of Brine Shrimp to 600ml tap water).
2. Using an airstone, bubble air into the water until all the salt has dissolved.
3. Add the brine shrimp eggs to the aerating solution. The eggs should hatch in 24-36 hours at a temperature of 27°C, or in 48 hours at 21°C.
4. To remove the live shrimp, remove the airstone and tilt the solution at an angle. The live shrimp will collect at the bottom of the container and the egg shells will float to the surface. Collect the live shrimp by siphoning into a very fine net.
5. Wash thoroughly in tap water before feeding to fish, axolotls, hydra, etc.

Ref: L5.00 20/03/09